

CoolCell™

Consistent -1°C/min cell freezing without alcohol

- Consistent -1°C/min cooling rate
- No alcohol
- Fast recycle time between freezes - minutes
- Unlimited number of freezing cycles - no maintenance
- No stuck lids, no frozen fingers
- Transportable without cold hands when frozen
- Unlike alcohol coolers, no warming of nearby freezer samples
- Eliminates the expense and disposal of isopropanol
- Holds up to 12 screw cap 2.0ml cryotube samples





The problem: Convenient cell freezing. Controlled -1°C/min cell freezing is currently a hassle. Alcohol-filled coolers need continuous maintenance, create hazardous waste, freeze fingers, don't open easily, and slow down work flow. No longer.

The solution: CoolCell. Achieve a consistent -1°C/min freeze rate without the use of hazardous solvent with CoolCell.

The solid state core and insulated design precisely balance heat removal during freezing to ensure repeatable, constant cooling all the way down to cryogenic storage temperature.

Cleaner, greener, easier to use.

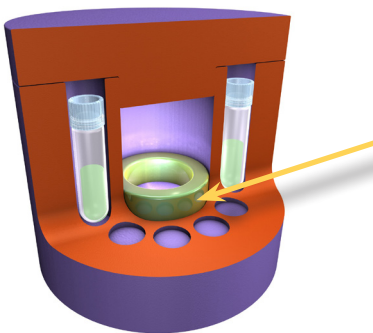

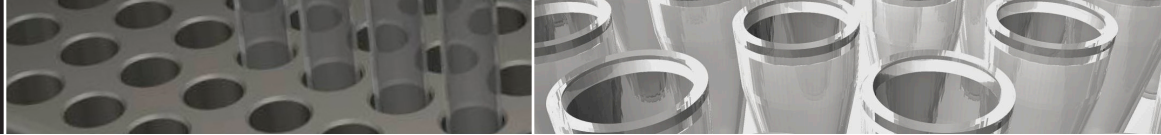

Save Cost, Stay Green: Factor the cost of 100% isopropanol, the time to decant and refill, the cost of hazardous material disposal, and the CoolCell makes immediate sense. Just say no to alcohol.



Always a constant -1°C/min: The CoolCell solid core design never needs changing and never varies in cooling rate. The insulated design provides repeatable, constant cooling throughout the freezing period. The result is a reliable -1°C/min rate every time.




Quick recycle time: Alcohol based freezers take a long, long time to warm up between cycles. Why wait? The CoolCell is ready to use “off the shelf” and is ready to use again just minutes after a freeze cycle. Eliminate waiting with the CoolCell.



Easy to handle: The insulated design protects your hands when you take the CoolCell out of the freezer. No more frosty hands. No more stuck lids.

Validation: The CoolCell and a standard alcohol filled freezer were compared to the ideal cooling rate of 1°C per minute (green).

Experiment: The temperature of a 1 ml volume of DMEM media, 10% DMSO, 20% fetal calf serum in a 2ml cryogenic storage vial was recorded using a thermocouple probe. Both the CoolCell and alcohol-filled freezing modules were filled to capacity with identical cryogenic vials and liquid loads. Freezing modules were placed into a -75°C freezer space. The plateau near 0°C reveals the latent heat released during the conversion of water to ice.

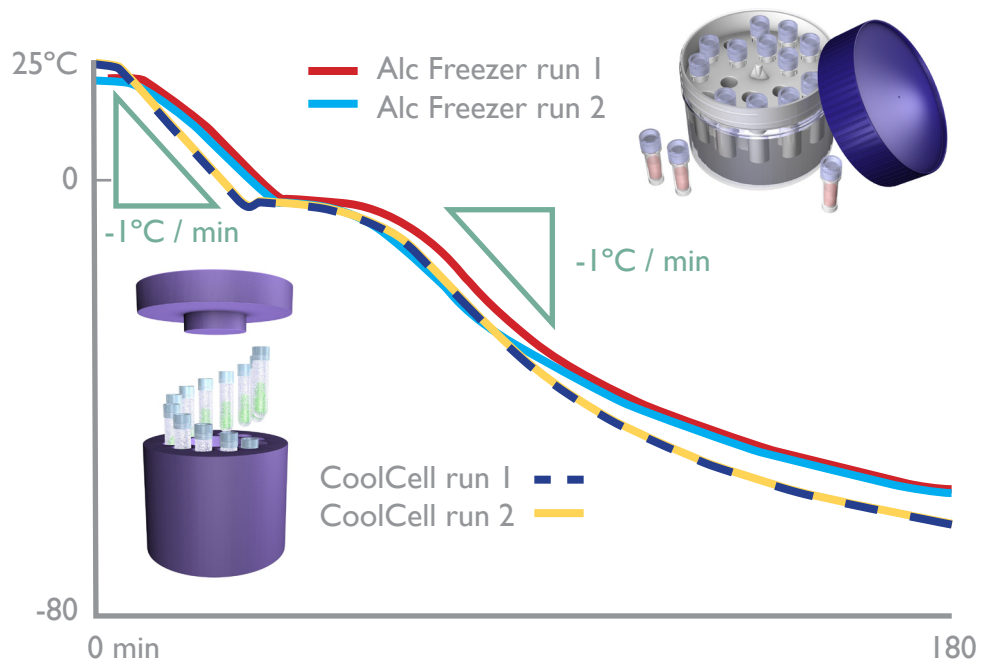

Conclusion: CoolCell provided superior freezing rate control when compared to an alcohol-filled freezing container and provided identical freezing profiles for each vial.

BCS-136

CoolCell

-1°C/min controlled rate freezing module for 1ml and 2ml cryotubes.
Capacity 12 tubes. Dimensions: 11Hx11.6dia cm Weight: 450g

Available from

UK
9 The Office Village
North Road, Loughborough
Leicestershire LE11 1QJ
United Kingdom
Tel. +44(0)1509 265265
Fax. +44(0)1509 269770
Email: sanyo@sanyo-biomedical.co.uk
www: www.sanyo-biomedical.co.uk

Netherlands
Nijverheidsweg 120
4879 AZ Etten Leur
The Netherlands
Tel: +31 (0)76 543 38 33
Fax: +31 (0)76 541 37 32
Email: welcome@sanyo-biomedical.nl
www: www.sanyo-biomedical.com

France Office
44, avenue de Valvins, BP 44
F-77212 Avon Cedex
France
Tel. +33 1 60719911
Fax. +33 1 60711693
Email: sanyo@wanadoo.fr



SANYO E&E Europe BV
Medical Division
Biomedical Business Unit